

## SERVICE NOTE

### JR6: Automatic holder adjustment

#### Procedure

- To adjust proper rotation of automatic holder, prepare screwdriver which is used for fixing of samples.
- Remove the automatic holder from the coils. Then turn with a screw to set proper mechanical tension of the inner shell (Fig. 1).
- The inner spherical shell of the holder should rotate freely but it should not be loosen too much.
- If you need to clean the automatic holder, unscrew the screw totally and pull out all other parts, as you can see in Fig. 2.
- We recommend to rinse the holder in water with small amount of detergent. Be careful to dry up all parts of the holder. **Using of alcohol to any part of holder is strictly forbidden.**



Figure 1: Adjustment of automatic holder



Figure 2: Adjusting screw