

SERVICE NOTE

JR: How to insert a holder

In the normal operation, the mechanical part of spinner magnetometer JR-6 is not visible. Installation of a holder to main shaft must be done without visual contact with main shaft. Following pictures illustrate how to insert a holder into the magnetometer:


1. This is a picture of main shaft with pin for leading of the holder:



2. Turn the holder until the drive pin enter the notch of the holder:



3. Continue pressing the holder downward until you overcome the pressure of an elastic pad under the shank of the holder.
4. Then continue clockwise turning in horizontal direction of bayonet groove until the end of horizontal part:

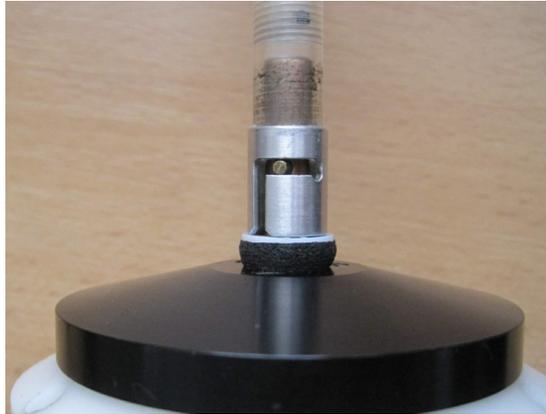

5. Finish the turning by upward jump of the holder which indicate proper installation of the holder:

